Mikroskop · Microscope ·

Duolux / BioDiscover / Biolux / Erudit / Biorit

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EN Instruction manual

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Contents

1 Imprint	12
2 Validity information	12
3 About this Instruction Manual	12
4 Intended Use	12
5 General safety instructions	13
6 Parts overview	15
7 Site location	16
8 Setting up power supply and turning on the device	16
9 Inserting slides	16
10 Observation.....	16
11 Cleaning and maintenance	17
12 Disposal.....	17
13 Warranty	18
14 EC Declaration of Conformity	18

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<http://www.bresser.de>

For information regarding liability claims or service requests, please refer to chapters "Warranty" and "Service" in this documentation. We ask for your understanding, that any requests or submissions that have been addressed to the manufacturer's address can not be processed.

Errors reserved - technical specifications subject to change.

Copyright: 2017Bresser GmbH

All rights reserved.

It is not allowed to reproduce this documentation or parts of it in any kind (e.g. photocopying, printing, etc) without written a permission of the manufacturer. It is also prohibited to use or transmit this documentation with electronic systems (e.g. image file, website, etc).

The descriptions and brand names of the corresponding firms used in this documentation are generally protected under trademark, brand and/or patent law for Germany, for the European Union and/or other countries.

2 Validity information

This documentation is valid for the products with the article numbers listed below:

5116200 8855000 8855001 8855002 8855003 8855004 8855100 9039000 9039001 5012000
5013000 5110000 5101000 micro-bio-beginner

Manual version: v062017a

Manual description: Manual_Biological-Microscopes-Beginners_de-en_BRESSER_v062017a

With any service inquiries, please state these information.

3 About this Instruction Manual

NOTICE

These operating instructions are to be considered a component of the device.

Please read the safety instructions and the operating instructions carefully before use.

Keep these instructions for renewed use at a later date. When the device is sold or given to someone else, the instruction manual must be provided to the new owner/user of the product.

4 Intended Use

- This device is not intended for use by individuals (including children) with limited physical, sensory or mental capabilities or those lacking in experience and/or knowledge, unless they are supervised by an individual responsible for their safety or have received instructions from them regarding the use of the device.
- This device is intended only for private use.
- It was developed for the magnified display of things in nature.
- The device is intended only for indoor use.

5 General safety instructions

DANGER

Danger of an electric shock!

This device contains electronic components that operate via a power source (power supply and/or batteries). In case of any improper use of this device, there is a risk of an electric shock. An electric shock can cause severe injury or even death. Therefore please read the safety instructions below to avoid an explosion.

- Disconnect the device from the power supply by pulling the power plug when it is not used or in case of longer interruption of operation and before starting any work on maintenance and cleaning.
- Position your device so that it can be disconnected from the power supply at any time. The wall socket you use should be located near the device and easily accessible, since the plug on the power cable serves as a disconnecting device for the power supply.
- Always pull on the plug to separate the device from the power supply. Never pull on the cable.
- Before operating, check the device, cables and connections for damage.
- Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.
- Only use the device in complete dry environment and do not touch it with wet or moist parts of your body.

DANGER

Danger of choking!

In case of any improper use of this device, there is a risk of choking, especially for children. Therefore please read the safety instructions below.

- Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard!
- This product contains small parts that could be swallowed by children. There is a risk of choking!
- If small parts has been swallowed, consult a doctor immediately!

DANGER

Danger of explosion!

In case of any improper use of this device, there is a risk of an explosion. Therefore please read the safety instructions below to avoid an explosion.

- Do not expose the device to high temperatures. Use only the supplied power adapter. Do not short-circuit the device or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.

CAUTION

Danger of injury!

This device contains components and/or accessories that can cause minor to severe injuries in case of any improper use. Therefore please read the safety instructions below to avoid any bodily injury.

- Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.

-
- Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after using the chemicals. In the event that the chemicals come into contact with your eyes or mouth, rinse thoroughly with water. If you are in pain after exposure, contact a doctor immediately and take the substances with which you came into contact with you.

CAUTION

Fire hazard!

In case of any improper use of this device, there is a risk of fire. Therefore please read the safety instructions below to avoid the initiation of burning.

-
- Never cover the ventilation slots or cooling fins of the device while using it or as long as it has not sufficiently cooled down!

NOTICE

Risk of property damage!

In case of any improper use of this device and/or its accessories, there is a risk of property damage. Therefore only use the device according to the safety instructions below.

-
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
 - Do not expose this device to higher temperatures and protect it from water and high humidity.
 - Protect the device from severe shocks!
 - For this device only use accessories and spare parts that comply with the technical information.

6 Parts overview

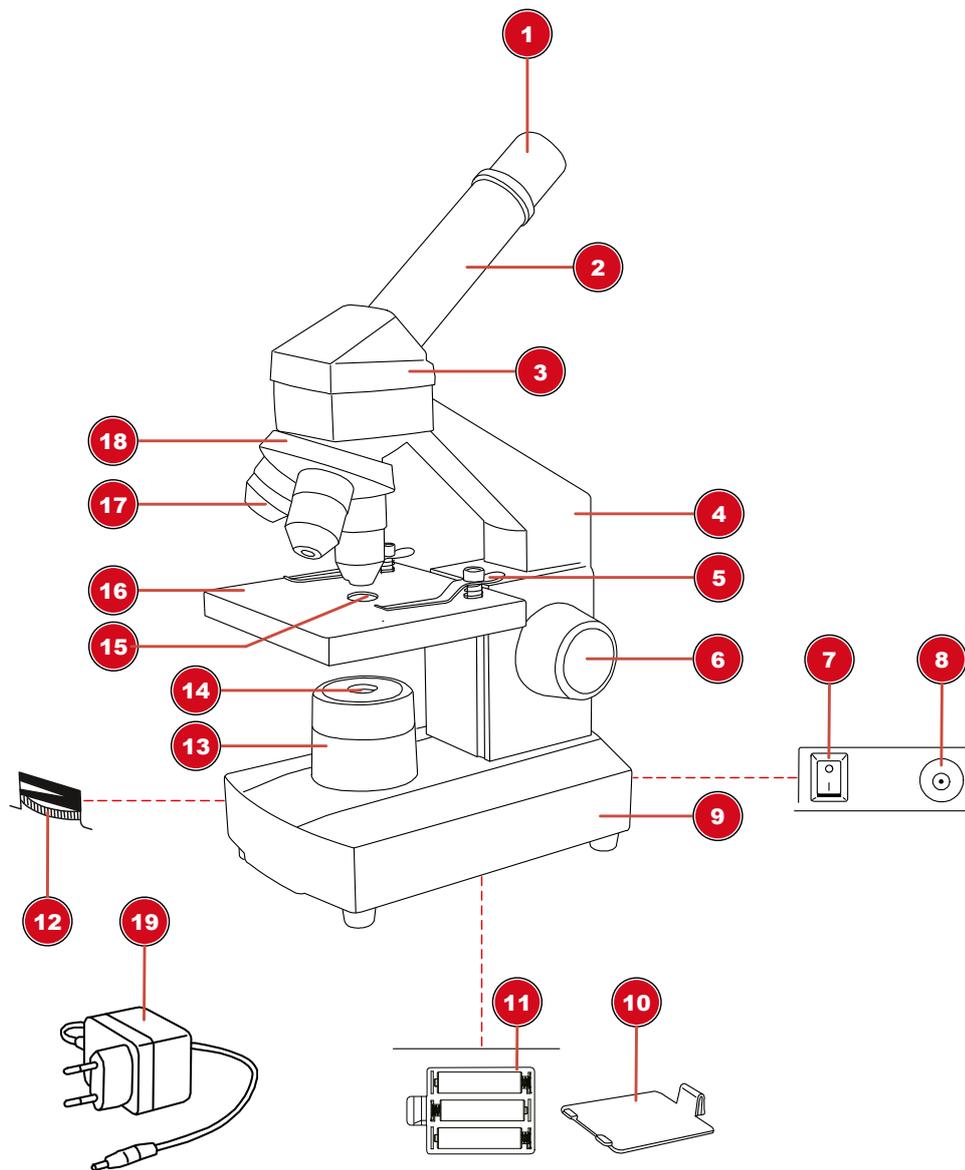
Illustration 2: Parts overview for biological microscopes

1 Eyepiece (can also be equipped with a pair of eyepieces) (optional equipment, depending on model)	2 Eyepiece holder
3 Microscope head	4 Microscope arm
5 Object fixing clamps	6 Focussing wheel
7 On/Off switch	8 DC socket for coaxial/barrel connector (can also be equipped with a C14 mains plug alternatively) (optional equipment, depending on model)
9 Microscope base	10 Battery compartment cover (optional equipment, depending on model)
11 Battery compartment (optional equipment, depending on model)	12 Rotary dimmer (optional equipment, depending on model)
13 Lighting unit	14 Lens
15 Light opening	16 Object stage
17 Objective	18 Objective nosepiece

19 Power adapter with Euro-plug, power cable and coaxial/barrel connector (can also be equipped with a Euro-plug and C13 mains socket) (optional equipment, depending on model)

7 Site location

Choose a suitable location before assembly and commissioning.
Place the device on a stable and flat surface, free from vibrations.

8 Setting up power supply and turning on the device

Depending on the model, the power supply can be obtained by battery or mains power (power adapter/cable).

Battery operation (depending on model)

1. Remove the battery compartment cover.
2. Insert the batteries into the battery compartment.
3. Re-attach the battery compartment cover.
4. Move ON/OFF switch to position 'I' to turn on the device.
5. Turn the rotary dimmer to set the illumination brightness. (optional function, depending on model)
6. Move On/Off switch to 'O' position to turn off the device.

Mains operation (depending on model)

1. Put the DC connector of the power cable into the DC connection socket of the device.
2. Put the Euro-plug of the power cable into the wall socket (230V maximum voltage).
3. Move ON/OFF switch to position 'I' to turn on the device.
4. Turn the rotary dimmer to set the illumination brightness. (optional function, depending on model)
5. Move On/Off switch to 'O' position to turn off the device.

9 Inserting slides

1. Move the object plate downwards completely by turning the focus wheel.
2. Push down both object clamps at the backside that they lift up at the frontside.
3. Slide the microscope slide from the frontside below the raised object clamps.
4. Let the object clamps fall back carefully until the slide is fixed under them.

10 Observation

NOTICE

Using eyepieces and objectives

When using a specific eyepiece or objective, you will get the individual magnification of an object being observed. The magnification is imprinted on the eyepiece and objective with a large number (e.g. 10x for 10-fold magnification, 4 for 4-fold magnification).

1. Move the object plate downwards completely by turning the focus wheel.
2. Insert a slide and place it in the middle under the objective.

3. Set an objective with lowest magnification below the object plate.
4. Insert an eyepiece with lowest magnification in the eyepiece holder.
5. Turn on the device and the illumination.
6. Look through the eyepiece and move the object plate upwards by turning the focus wheel until a sharp image is visible.
7. If necessary, adjust the image sharpness by turning the fine focus adjustment.

NOTICE

Avoid damages to the specimen!

Make sure you turn the focus wheel very slowly, to avoid the objective touch the slide or specimen, and thus to avoid damages to either of them, or to both. Only turn the focus wheel until it stops and do not apply excessive pressure.

11 Cleaning and maintenance

- Protect the device against dust and moisture.
- Avoid putting fingerprints and equal contaminations on any optical surfaces.
- Make sure the residual moisture is fully eliminated before use.
- Disconnect the device from the power supply by pulling the power plug when it is not used or in case of longer interruption of operation and before starting any work on maintenance and cleaning.
- Only use a dry cloth to clean the exterior of the device. To avoid damaging the electronics, do not use any cleaning fluid.
- Do not use organic solvents (e.g. alcohol, ether, acetone, xylol or other dilutions) to clean lacquered parts or plastic parts!
- Disassemble optical elements (e.g. objective, eyepiece, etc) before cleaning.
- Blow away loose dust from the lens surfaces first.
- Use high quality lens cleaning towels or soft cloth and moisten it with a little bit of pure alcohol (available in a drug store). Wipe the lens surface with it.
- Pull a dust protection cover* over the microscope and store it in a closed container at a dry and mould-free place. Store the microscope and the accessories in the relevant containers when they are not used for a longer time. It is recommended to store also objectives and eyepieces in closed containers with desiccant.

(*depending on the model a dust protection cover and box can be part of the content)

NOTICE

Remember:

A well maintained microscope will keep its optical quality for years and thus maintain its value.

12 Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please make sure to dispose of your used batteries as required by law — at a local collection point or in the retail market. Disposal in domestic waste violates the Battery Directive.

Batteries that contain toxins are marked with a sign and a chemical symbol. "Cd" = cadmium, "Hg" = mercury, "Pb" = lead.

13 Warranty

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at www.bresser.de/warranty_terms.

14 EC Declaration of Conformity

 Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards.

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GB IE

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

e-mail: service@bresseruk.com
Telephone*: +44 1342 837 098

BRESSER UK Ltd
Unit 1 starborough Farm,
Starborough Road, Nr Marsh Green,
Edenbridge, Kent TN8 5RB
Great Britain

*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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Folux B.V.