# Levenhuk Kelvin HPM25 Handheld Photomicroscope

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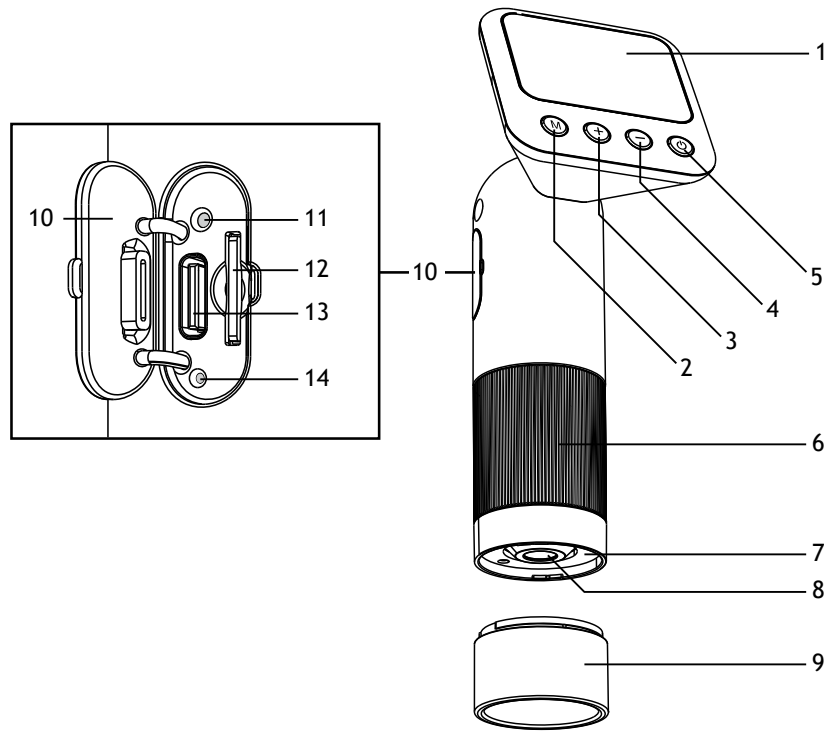
# EN Levenhuk Kelvin HPM25 Handheld Photomicroscope

**Caution! Never look directly at the Sun through this device, as this may cause permanent eye damage and even blindness.**

The kit includes: digital photomicroscope, USB Type-C cable, pouch, cleaning cloth, user manual, and warranty.

## Charging the device

This device uses a rechargeable Li-ion battery. Connect the USB Type-C cable to the USB Type-C slot (13) and power bank, PC, or power adapter (not included). The power indicator (11) will light red. Charging takes approximately 2 hours. When charging is complete, the power indicator will light green.

## Getting started

- Press and hold (5) to turn the device on. The LED illumination (7) will light up.
- Press and hold (5) to turn the device off.

## Installing/removing the memory card

- Insert the microSD card (not included) into the memory card slot (12). Make sure the direction is right. Do not force the card in.
- Format the microSD card before starting, as it will improve its compatibility with the device.
- To remove the card from the slot, press the edge of the memory card in gently, and the card will pop out. Then, remove the card.

## Mode selection

Press (2) to change the mode. The setting order: Photo mode > Video mode > Playback mode. The default mode is photo mode.

## Settings

Press and hold (2) to enter settings, then press (2) to choose the settings menu. The setting order: Photo settings > Video settings > General settings. Use (3) to scroll up and (4) to scroll down, then press (5) to confirm or (2) to return. You may now use the device to observe objects, take photos, and record videos.

## Usage

- Place the object to be observed on a flat surface.
- Ensure that the object is correctly positioned inside the transparent plastic tube (9) below the objective lens (8) (Fig. 1).

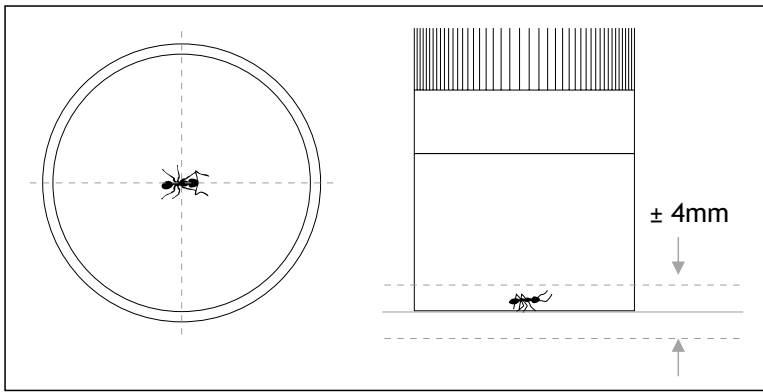
Fig. 1

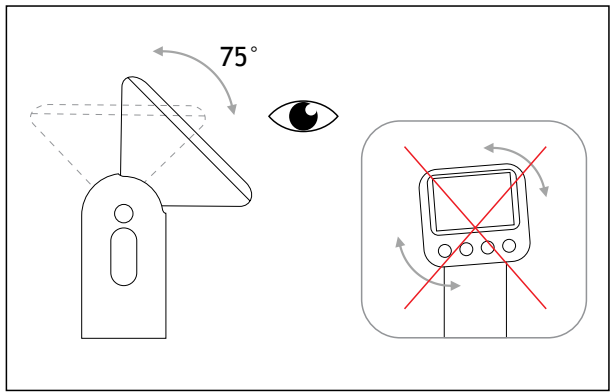
Fig. 2

- Tilt the screen (1) to a comfortable viewing angle. Please note that the screen is not to be rotated horizontally (Fig. 2).
- Rotate the focusing ring (6) until the view is in focus.

## Photo and video

Press (5) to take a photo or to start recording a video. Press (3) to zoom in and (4) to zoom out. Press and hold (3) to turn the LED illumination (7) on and increase brightness, and (4) to decrease brightness and turn the LED illumination off.

## Playback

Press (3) or (4) for scrolling through your pictures and video. Press (5) to playback a video file, then press and hold (3) or (4) to fast forward or rewind the video. Press and hold (2) to delete the current file.

## Copying files

You can transfer files from the device to your PC directly.

- Insert the memory card into a card reader; or
- Connect the device to your computer with a USB Type-C cable.

## Software

To work with the device on a PC, you need to install the VideoCap software, available for download on the official Levenhuk website. This software enables you to view your photos and videos in real time on your PC screen.

## Specifications

Research method	bright field
Magnification	400–800x
Min. size of the objects to be seen	3µm
Field of view	2.3mm
Optics material	optical glass
Optics coating	multi-coated
Objective lens type	achromatic
Microscope head	rotatable digital screen (up to 75°) / PC screen
Display	2" color IPS screen, 320x240px
Sensor	CMOS, 720p
Photo format and resolution	*.jpg: 1MP, 2MP
Video format and resolution	*.avi: 720p, 1080p
Brightness level	3 modes (high/low/off)
LED illumination	6 pcs white LEDs
Display auto-off	30s, 1 min, 2 min
Device auto-off	3 min, 5 min, 10 min
Storage	microSD card up to 128GB (not included)
Software requirements	Windows 7 and above
Operating temperature range	-10... +40°C (+14... 104°F)
Power supply	3.7V, 700mA·h rechargeable Li-ion battery; 5V, 1A DC adapter (not included); USB Type-C cable (included)
Battery duration	2–2.5h

The manufacturer reserves the right to make changes to the product range and specifications without prior notice.

## Care and maintenance

Take the necessary precautions when using the device with children or others who have not read or who do not fully understand these instructions. Do not try to disassemble the device on your own for any reason. Protect the device from water, dust, sudden impact and excessive mechanical force. Do not touch the optical surfaces with your fingers. Clean the lens surface with compressed air or a soft lens cleaning cloth. Never attempt to operate a damaged device or a device with damaged electrical parts! Store the device in a dry cool place. **If a part of the device or battery is swallowed, seek medical attention immediately.**

## Battery safety instructions

The device is equipped with a rechargeable lithium-ion battery. This avoids frequent battery replacement. Always switch the device off when not in use. If the battery charge is low, please recharge the device in time. Do not overheat the battery. Do not discharge the battery completely. Keep batteries out of the reach of children, to avoid risk of ingestion, suffocation, or poisoning. Utilize used batteries as prescribed by your country's laws.

## Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars, and other optical products, except for their accessories, carry a **lifetime warranty** against defects in materials and workmanship. A lifetime warranty is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for **six months** from the purchase date. The warranty entitles you to the free repair or replacement of the Levenhuk product in any country where a Levenhuk office is located if all the warranty conditions are met.

For further details, please visit: [levenhuk.com/warranty](http://levenhuk.com/warranty)

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch.

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Levenhuk® is registered trademark of Levenhuk, Inc.
© 2006–2026 Levenhuk, Inc. All rights reserved.
Levenhuk Inc. (USA): 6021 Catlin Dr., Tampa, FL 33612, USA,