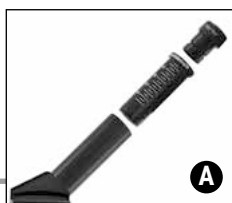
MIKROSKOP MICROSCOPE

40X-1024X



- DE** Schnellstart-Anleitung
- EN** Quick-start instructions
- FR** Guide de démarrage rapide
- NL** Snelstart-handleiding
- IT** Guida rapida
- ES** Guía de inicio rápido
- RU** Краткое руководство пользователя


Lieferumfang · Contents · Contenu de la livraison · Inhoud · Accessori in dotazione
Artículos incluidos en la entrega · Комплект поставки

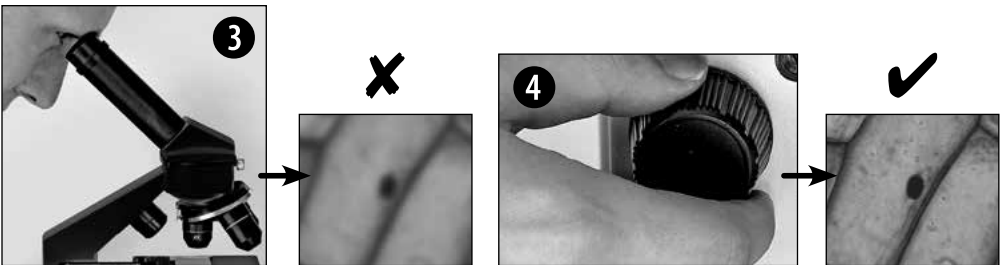
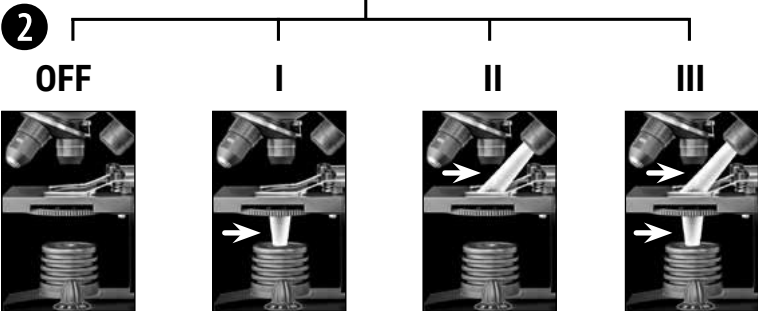
Stromversorgung / Beleuchtung · Power supply and illumination operation · Alimentation électrique / Eclairage · Stroomvoorziening / Belichting · Alimentazione elettrica / Illuminazione · Alimentación de corriente / Iluminación · Питание от сети / Подсветка

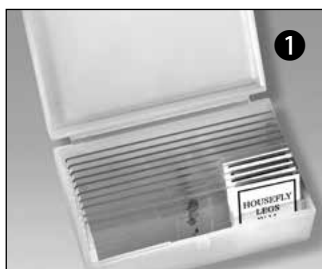
Zubehör anbringen · Accessory installation · Fixation des accessoires · Accessoires aansluiten ·
Montaggio degli accessori · Instalación de accesorios · Установка дополнительного оборудования

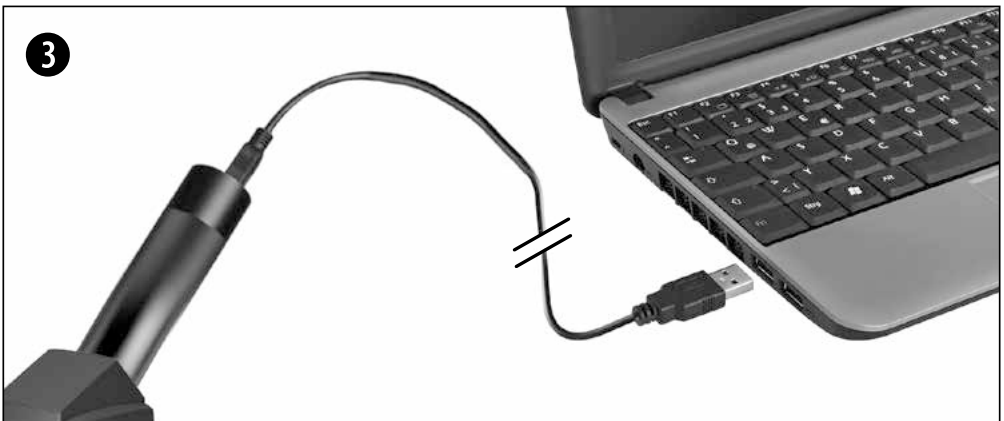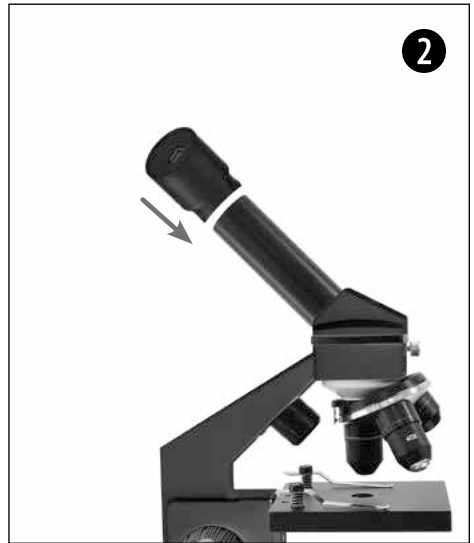PC-Okular anschließen · Installing the PC eyepiece · Raccordement de l'oculaire PC · PC-oculair aansluiten · Collegamento dell'oculare PC · Conexión del ocular para PC · Установка PC окуляра




4





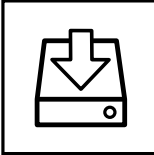
SOFTWARE DOWNLOAD:

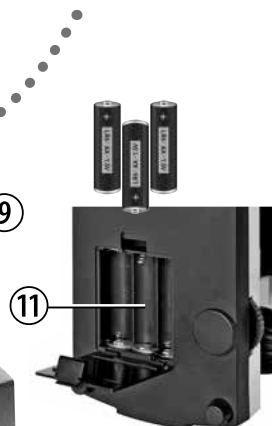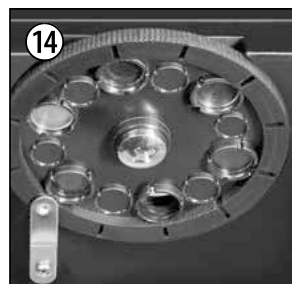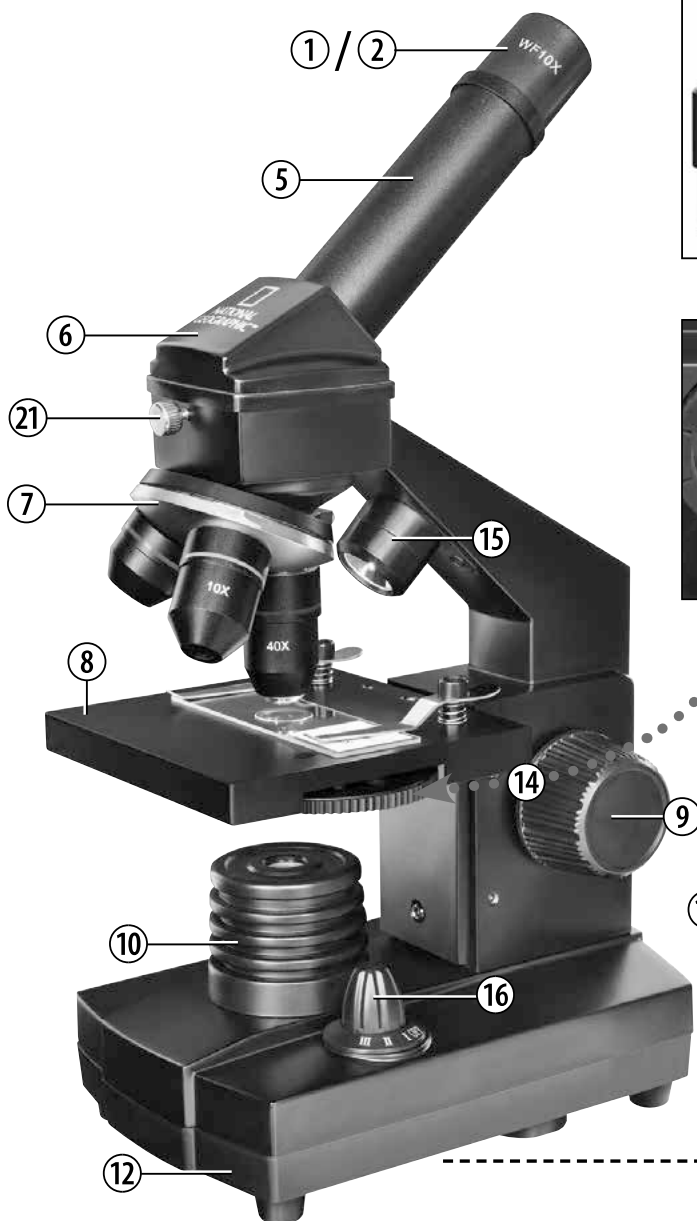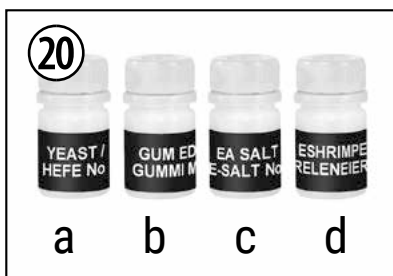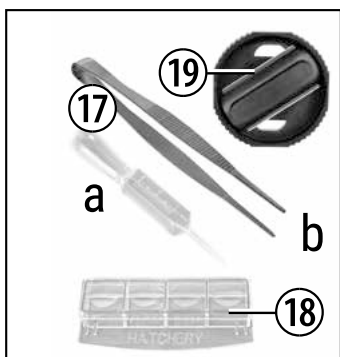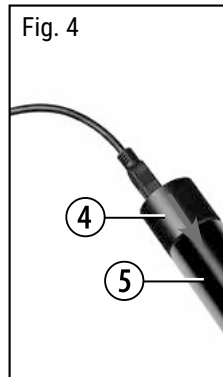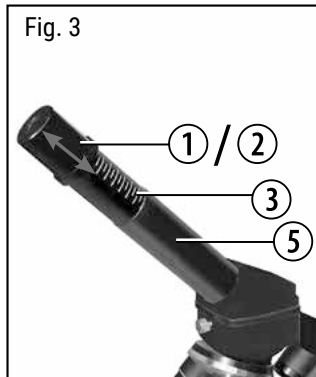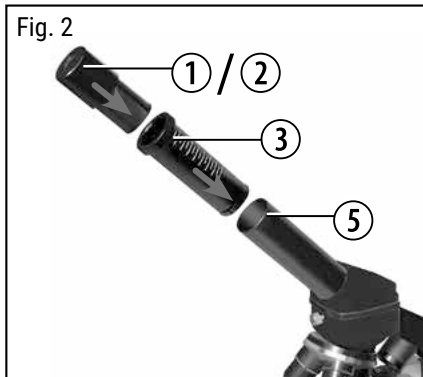www.bresser.de/download/CamLabLite

Download: <http://www.bresser.de/9039100>

Eine ausführliche Anleitung steht für Sie auf unserer Internetseite zu Ihrer Verfügung.
A more detailed manual is available on our website. • Un guide détaillé est disponible sur notre site. • Een uitgebreide handleiding is beschikbaar op onze website. • Un manuale di istruzioni dettagliato è disponibile sul nostro sito Web su internet. • Un manual de Instrucciones detallado se encuentra disponible en nuestra página Web en internet.
Подробное руководство можно найти на веб-сайте.

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EN About this manual

This quick-start guide will help you put the device into operation quickly, but it is not intended to replace the more detailed user manual.

Please read the safety instructions and the user manual carefully before using the device. This quick-start manual contains the most important steps to start using your new device. Comprehensive operating instructions and further information, including troubleshooting advice, is available on our website: www.bresser.de (Download: <http://www.bresser.de/9039100>)

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General Warnings

- **Choking hazard** – This product contains small parts that could be swallowed by children. This poses a choking hazard.
- **Risk of electric shock** – This device contains electronic components that operate via a power source (batteries). Only use the device as described in the manual, otherwise you run the risk of an electric shock.
- **Risk of fire/explosion** – Do not expose the device to high temperatures. Use only the recommended batteries. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.
- **Risk of chemical burn** – Make sure you insert the batteries correctly. Empty or damaged batteries could cause burns if they come into contact with the skin. If necessary, wear adequate gloves for protection.
- Leaking battery acid can lead to chemical burns. Avoid contact of battery acid with skin, eyes and mucous membranes. In the event of contact, rinse the affected region immediately with a plenty of water and seek medical attention.
- Use only the recommended batteries. Always replace weak or empty batteries with a new, complete set of batteries at full capacity. Do not use batteries from different brands, types or with different capacities. Remove the batteries from the unit if it is not to be used for a long period of time, or if the batteries are empty!
- Never recharge normal, non-rechargeable batteries. This could lead to explosion during the charging process.
- Rechargeable batteries are only to be charged under adult supervision.
- Rechargeable batteries are to be removed from the toy before being charged
- The terminals must not be short-circuited.
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Tools with sharp edges are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.
- Keep instructions and packaging as they contain important information.

EC Declaration of Conformity

Bresser GmbH has issued a „Declaration of Conformity“ in accordance with applicable guidelines and corresponding standards. The full text of the EU declaration of conformity is available at the following internet address:

www.bresser.de/download/9039100/CE/9039100_CE.pdf

Parts overview:

- ❶ 10X WF eyepiece
- ❷ 16X WF eyepiece
- ❸ Barlow lens
- ❹ Electronic eyepiece (MicrOcular)
- ❺ Microscope head
- ❻ Objective revolver
- ❼ Microscope stage
- ❽ Focus wheel
- ❾ LED lighting (transmitted light)
- ❿ Electricity supply
- ⓫ Microscope base
- ⓬ Main plug
- ⓭ 5 slides, 10 covering glasses and 5 preparations in a plastic box
- ⓮ Colour filter disc

- 15 LED lighting (reflected light)
- 16 Direct light/transmitted light switch
- 17 Microscope tools: a) pipette; b) tweezers
- 18 Hatchery
- 19 MicroCut
- 20 Specimens: a) yeast; b) gum media (specimen inclusion medium); c) sea salt; d) brine shrimp eggs
- 21 Locking screw

How do I install the MicrOcular?

Note:

The MicrOcular only works without the included Barlow lens! The magnification setting is changed when the MicrOcular is used, and it must be readjusted with the focus knob.

First, remove the Barlow lens with the current eyepiece from the eyepiece support and place the MicrOcular in the eyepiece.

How do I install the software?

1. You can download the setup file for installation here:
www.bresser.de/download/CamLabLite
2. Start the installation with a double click on the .exe file and follow the instructions of the installation manager.
3. It is not necessary to install a special driver. The camera is automatically recognized by your Windows operating system.

Using the MicrOcular

Preparation

1. Slide a specimen under your microscope and focus on it.
2. Remove the eyepiece and Barlow lens from the eyepiece support and the dust cap from your MikrOkular and then install the MikrOkular into the eyepiece support instead of the Barlow lens.
3. Start your PC if you haven't yet and connect your MikrOkular to the USB port of your computer.
4. Start the „CamLabLite“ software and activate the camera by clicking on „USB Video Device“.
5. Click on „Snapshot“ to take a picture or on „Capture“ to record a video.
6. Click on the floppy disk icon to save the image or capture.

Experiments

Use the following web link to find interesting experiments you can try out.

<http://www.bresser.de/downloads>

Warranty & Service

The regular guarantee period is 5 years and begins on the day of purchase. You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at www.bresser.de/warranty_terms.

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© National Geographic Partners LLC. All rights reserved.
trademarks of National Geographic Society, used under license.