LEVENHUK ERGO 40 PRO DIGITAL MAGNIFIER

EN User Manual

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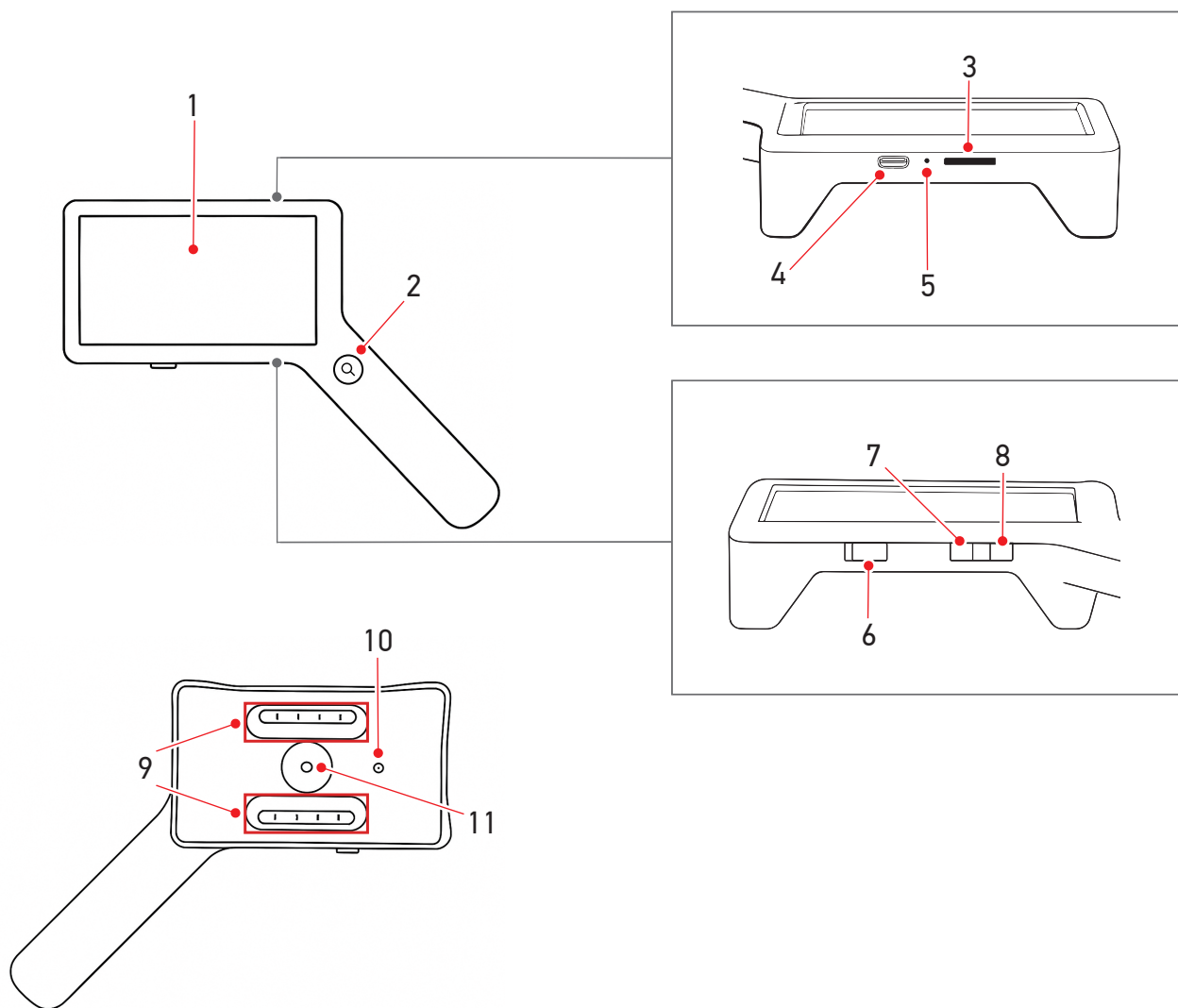
EN Levenhuk ERGO 40 PRO Digital Magnifier

Never, under any circumstances, look directly at the Sun, another bright source of light or at a laser through this device, as this may cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS.

The kit includes: digital magnifier, USB Type-C cable, user manual, and warranty.

Charging the device

This device uses a rechargeable Li-ion battery. Connect the USB Type-C cable to the USB Type-C slot (4) and power adapter (not included) or PC. The power indicator (5) will light red. Charging takes approximately 2 hours. When charging is complete, the power indicator will light green.

! If the low battery indicator appears on the screen (1), charge the device.

Getting started

- Turn the power switch (6) on. The white LEDs (9) will light up. Place the object to be observed directly below the objective lens (11).
- Turn the power switch (6) off. The device will turn off automatically if it is not in use for 10 minutes. Press (2) to turn it on again.
- If the screen freezes or the device fails to start, it must be charged or restarted. Press (10) using a thin rod to restart the device. Then press (6) to turn it on again.

Installing/removing the memory card

- Insert the microSD card (not included) into the microSD card slot (3). Make sure the direction is right. Do not force the card in.
- Format the microSD card before starting, as it will improve its compatibility with the device.
- To remove the card from the slot, press the edge of the memory card in gently, and the card will pop out. Then, remove the card.

Time setting

- Connect the device to a PC using a USB Type-C cable. Plug the USB Type-C connector of the cable into the USB Type-C slot (4), then plug the standard USB connector of the cable into the PC.
- Open the time.txt file on the USB drive and enter the current date and time. Date format: YYYY.MM.DD, time format: hh:mm:ss.
- Save the file, and the date and time will be displayed on the device.

Usage

Magnification setting

Press (2) to change the magnification. The setting order: 6x > 10x > 13x > 17x.

LED brightness setting

Press (7) to change the LED brightness. The setting order: 100% > 75% > 50% > Off.

Overlay setting

Press and hold (2) to show the overlay. The setting order: Reticle > Grid > Off.

Capturing photo

Press (8) to capture a photo. Keep the device still about 1 second after capturing. The captured photos will be automatically saved to the memory card. Press (8) again to return to normal mode.

Taking video

Press and hold (8) for 3 seconds to enter the video mode. Press and hold (8) for 3 seconds again to exit. The video will be automatically saved to the memory card.

Copying files

- Connect the device to a PC using a USB Type-C cable. Turn the power switch (6) on.
- Open the DCIM folder on the USB drive and navigate to the DCIMA folder to view the photo and video files. You can upload them to your PC.
- Turn the power switch (6) off.

Specifications

Display	4" color LCD, 800x480px
Sensor	1/4" CMOS, 1280x720px
Focus distance	15–19mm
Field of view	45°
Magnification	6x, 10x, 13x, 17x

LED illumination	4 pcs white LEDs
Storage	microSD card up to 128GB (not included)
Photo format and resolution	*.jpg, 1920x1080px
Video format and resolution	*.avi, 1280x720px
Auto-off	10 min
Operating temperature range	0... +40°C (+32... 104°F)
Power supply	3.7V, 1200mA-h rechargeable Li-ion battery; 5V, 1A DC adapter (not included); Type-C USB cable (included)
Battery duration	180 min

The manufacturer reserves the right to make changes to the product range and specifications without prior notice.

Care and maintenance

Take the necessary precautions when using the device with children or others who have not read or who do not fully understand these instructions. Do not try to disassemble the device on your own for any reason. Protect the device from water, dust, sudden impact and excessive mechanical force. Do not expose the device to direct sunlight or other light sources. Use only dry, lint-free cloths to wipe the outer parts of the device. Do not use cleaning fluid to avoid the electronics damage. To avoid scratching the lenses, only use a soft cloth and do not press too hard. Dry your hands thoroughly before using the device, otherwise it may result in electric shock. Never point the LED lights into your eyes. In case of malfunction, stop using the device. **For repairs and cleaning of any kind, please contact your local specialized service center.**

Battery safety instructions

The device is equipped with a rechargeable lithium-ion battery. This avoids frequent battery replacement. Always switch the device off when not in use. If the battery charge is low, please recharge the device in time. Do not overheat the battery. Do not discharge the battery completely. Keep batteries out of the reach of children, to avoid risk of ingestion, suffocation, or poisoning. Utilize used batteries as prescribed by your country's laws.

Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars, and other optical products, except for their accessories, carry a **lifetime warranty** against defects in materials and workmanship. A lifetime warranty is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for **six months** from the purchase date. The warranty entitles you to the free repair or replacement of the Levenhuk product in any country where a Levenhuk office is located if all the warranty conditions are met.

For further details, please visit: levenhuk.com/warranty

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch.

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Levenhuk® is a registered trademark of Levenhuk, Inc.
© 2006–2026 Levenhuk, Inc. All rights reserved.