S-Viewer Software

- EN User Manual
- BG Ръководство за потребителя
- CZ Návod k použití
- DE Bedienungsanleitung
- ES Guía de usuario
- HU Használati útmutató
- IT Guida all' utilizzo
- PL Instrukcja obsługi
- PT Manual do usuário
- RU Руководство пользователя

Levenhuk Inc. (USA): 6021 Cattin Dr., Tampa, FL 33612, USA,
+1 813 468-3001, contact_us@levenhuk.com
Levenhuk Optics s.r.o. (Europe): V Chotejně 700/7,
102 00 Prague 102, Czech Republic, +420 737-004-919, sales-info@levenhuk.cz
Levenhuk® is registered trademark of Levenhuk, Inc.
© 2006–2026 Levenhuk, Inc. All rights reserved.
levenhuk.com
20260626

levenhuk 

EN S-Viewer Software


S-Viewer is professional software for digital microscopes. It provides real-time video processing, dynamic measurements, object counting, image stitching, HDR, EDF, and flexible layer management. This guide covers the essential steps to get you started.

Running the application


1. Launch the software after installation. The first time you start the software, it will automatically detect a video capture device and show live preview.
2. If multiple cameras are connected, select the target device:
Menu → View → Camera control panel → choose the device from the list.
3. To change the language:
Menu → Option → Language → select desired language (restart required).

Main window interface

Interface elements: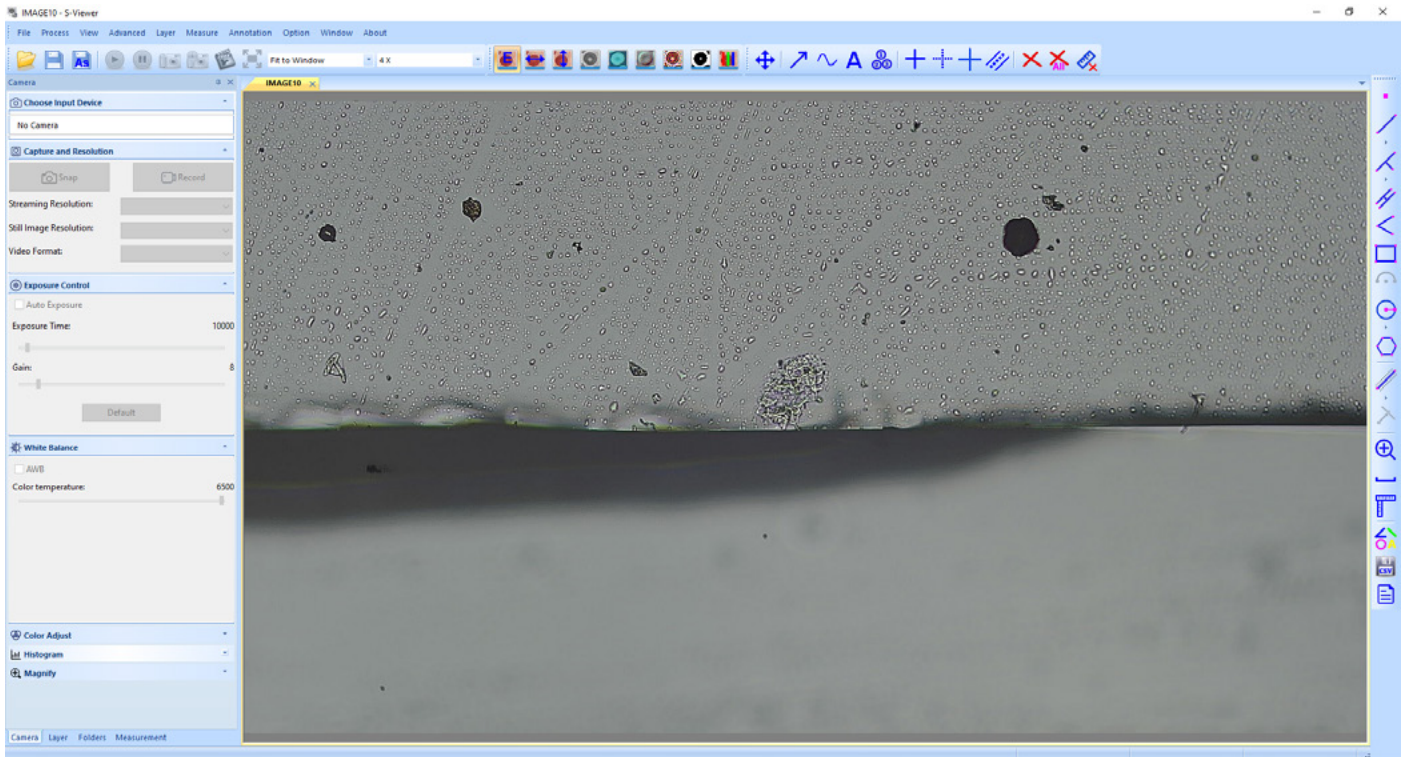


1. **Menu bar**
 Contains all the action commands for the application.

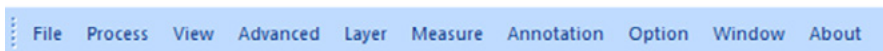2. **Toolbar**
 Contains buttons for the most commonly used actions.

3. **Toolbox tabs**
 Grants access to quick setup and editing tools.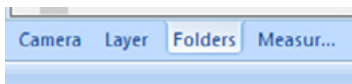

Name	Description
Camera	Select camera, capture/resolution, exposure, white balance, color adjustment, histogram, and magnifier
Layer	Manage layers (new, delete, rename, show/hide, set current)
Measurements	View geometry, coordinates, and appearance of measurement objects after moving them
Wi-Fi Camera	Search, connect and configure Wi-Fi microscope cameras (supports multiple simultaneous views)
Folders	Local file browser; double-click thumbnails to open images (BMP, JPG, PNG, TIF, GIF)


4. **Workspace**
 Contains all the opened images, ready for editing.

5. Processing toolbar

Contains tools and real time effects (flip, grayscale, emboss, sharpen, binary, negative, RGB match).

Camera menu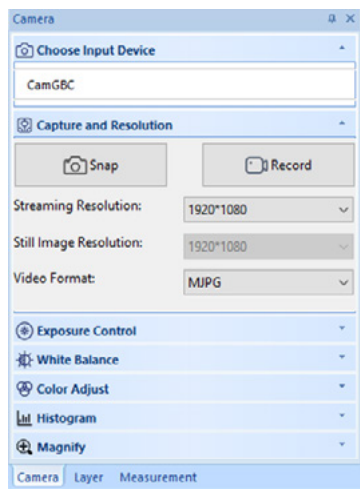


Viewing

Clicking on the camera name will allow you to see the live camera view on your PC monitor. A CamGBC tab will appear in the workspace.

You can edit the resulting image with Exposure Control, White Balance, Color Adjust and other tools on the Camera tab of the toolbox.

Capturing an image

Click the **Snap** button. The current view will be captured and the resulting image shown in a new workspace tab. Every time you click the Snap button, a new image is taken and shown in a new tab.

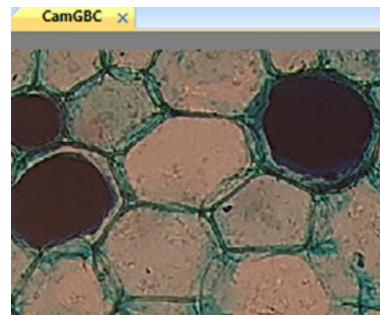Capturing a video

Click the **Record** button. Choose the file format, a name for your video and the output folder. Upon confirming these selections, the live feed from the camera will be saved as a video until you choose to stop it.

Editing the image

Saving

You can save one or several images on your hard drive by choosing the required option in the File menu. Choose between **Save** and **Save as...**

The following file formats are supported:

- Window Bitmap (*.bmp)
- JPEG (*.jpg)
- Portable Network Graphics (*.png)
- Tag Image File Format (*.tif)
- PCX (*.pcx)
- Targa (*.tga)

You can embed Title, Author, Date/Time, Ruler, and Measurement data into the image.

Measurements

S-Viewer supports dynamic (live video) and static image measurement.

Calibrate first: **Measure** → **Calibration** sets magnification, resolution, ruler unit, and actual length.

Measurement tools are in the **Measure** menu or the **Measurement** control panel. Measured data can be exported as CSV file.

Tool	Description
Point	Places a point and gives you its X and Y coordinates
Line → Draw Line	Draws a line between two points on a given layer
Line → Draw Horizontal Line	Draws a horizontal line between two points on a given layer
Line → Draw Vertical Line	Draws a vertical line between two points on a given layer
Line → Draw Polyline	Draws a polyline between many points on a given layer
Vertical → Draw Vertline	Draws a line perpendicular to the given line
Vertical → Draw Vertlines	Draws several lines perpendicular to the given line
Parallel	Draws two parallel lines and numbers them
Angle	Draws an angle and measures it
Draw Rectangle	Draws a rectangle from two points along a diagonal and shows its dimensions
Draw Arc	Draws an arc along three points and shows its length
Circle	Draws a circle on a given layer, using the Two Points method
Tri-circle	Draws a circle on a given layer, using the Three Points method
Circle to vertical line	Draws a circle on a given layer, using the Center + Radius method
Annulus	Draws a ring (two circles with a common center) and shows its inner and outer radius
Two Circles	Draws two circles (Three Points method) and a line connecting their centers. Shows the distance between these points
Polygon	Using your mouse, place as many points as you need. Right-click to create a polygon
Scale bar	Sets the scale according to the length and unit after calibration
Calibration	The default is 15 sets of magnifications, all of which need to be re-calibrated and can be renamed in the management magnification
Auxiliary magnification	Zooms in on the current area based on the mouse position
Measurement settings	Allows to set the scale, font size and background color of the measurement

The measurements of all the shapes on the image may be seen in **View → Measurement Sheet** menu.

Measurement Sheet											
Index	Name	Length	Distance	Width	Height	Center	Diameter	Area	Angle	Start	End
1	A1								295.207	(226.800,120.00...	(257.800,66.600)
2	Pt1					(136.400,58.000)					
3	Pt2					(80.400,129.200)					
4	Vt11	93.392							117.553		
5	Vt12	42.293							51.721		
6	Vt13	12.750							166.399		
7	Cv1	84.257				(246.000,25.600)		12948.502			
8	Tc1	310.389 233.734	43.953				98.800 74.400	5773.237 4347....			

Image processing

Access via **Menu → Process**.

Icon	Function	Description
	Flip horizontal/vertical	Mirrors image in real time
	Gray	Converts to black and white
	Emboss	Adds emboss effect
	Sharpen	Enhances edges
	Binary	Adjusts threshold (1–255) for black/white display
	Negative image	Inverts colors
	RGB adjustments	Adjusts Red/Green/Blue balance



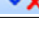
Layers

Default image opens as **background layer**.

Function	Description
New	Create named layer
Delete	Remove selected layer (cannot delete background or current layer)
Delete All	Keep only background
Current	Set active layer
Show/Hide	Hide any layer (background and current cannot be hidden)
Rename	Rename current layer

Annotations and overlays

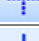Access via toolbar or **Menu → Annotation**.

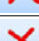Icon	Function	Description
	Select/Move	Moves added graphics/text
	Arrow	Draws an arrow
	Curve	Draws a freehand curve
	Text	Adds text
	Number markers	Adds serial numbers for counting
	X-Y guide solid	Adds solid crosshair lines
	X-Y guide dashed	Adds dashed crosshair lines
	Graduated scale	Adds a graduated scale
	Center line	Draws a center line between two parallel lines
	Delete	Removes an annotation
	Delete all	Removes all annotations
	No action	Cancel current measurement state

Mark overlay (via **Menu → Annotation**) opens overlay settings: choose circle/rectangle, set width/height/radius (as percentage of video window), X/Y offset.

Settings

To customize the application, you can use the **Options** menu.

Save Option: sets filename prefix, start suffix, folder path, image type (Bitmap/JPEG), and what to embed (Title, Date/Time, Measurement, Ruler, Description, Author).

UI Style: switches between different interface styles.

Language: changes application language (restart required).

Troubleshooting (previewing fails, blank screen)

1. Check the light source and exposure value.
2. Verify the video device connection and driver installation (**Device manager**).
3. Ensure the correct video device is selected (**View** → **Camera** control panel).
4. Close and restart the software.
5. If the problem persists, unplug and reconnect the camera, then restart the software.
6. If it does not help, restart your PC.

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