Photomicroscopes

Levenhuk Kelvin Touch

EN User Manual

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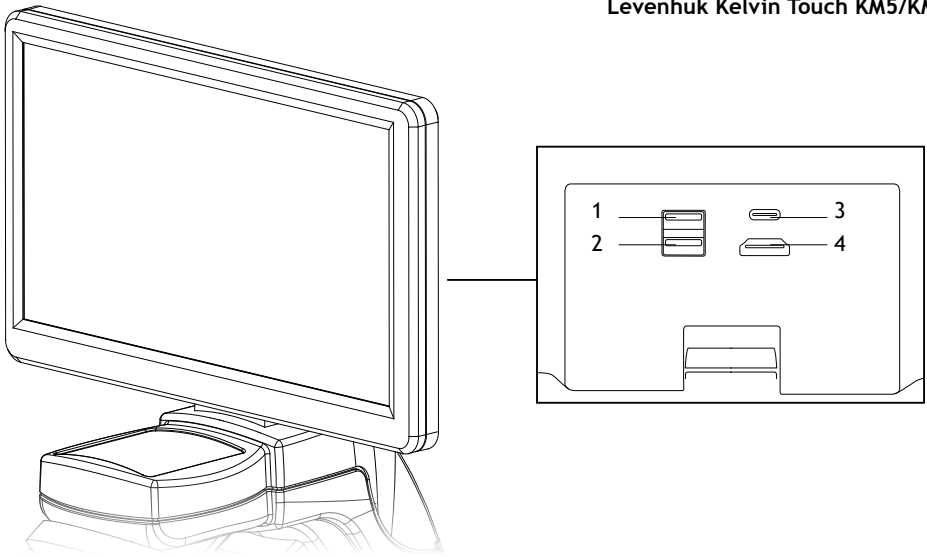
Caution! Please refer to the specifications table for the correct mains voltage and never attempt to plug a 110V device into 220V outlet and vice versa without using a converter. Remember that mains voltage in the U.S. and Canada is 110V and 220–240V in most European countries.

The kit includes: photomicroscope, power adapter, USB Type-C cable, microSD card, user manual, and warranty.

Getting started

- Plug the USB Type-C cable into the power cable socket on the microscope body (15) and connect it to a power source using the power adapter; or
- Open the battery compartment cover (16) and insert the battery (not included) according to the correct polarity.
- Press (14) to turn the device on/off.

Downloading the app (KM3 and KM7 models)

- Scan the QR code below to download the TSViewer app for iOS from the Apple App Store:

<https://apps.apple.com/us/app/nbtsview/id6448884238>

- Download the TSViewer app for Android from the official Levenhuk website.
- Enable the permissions for the app.

Pairing with a smartphone (KM3 and KM7 models)

- Press (14) to turn the microscope on. Open the Wi-Fi settings on your device and connect to the “levenhuk-XXX” Wi-Fi network.
- Start the app on your device, select the “levenhuk-XXX” Wi-Fi network, and tap “Connect”.

Usage

- Turn on the power and the illumination will turn on. Set the brightness at approximately 70%. You can use upper or lower illumination. Place the specimen on the stage. Make sure the clips hold the specimen firmly in place.
- To change the objective magnification, turn the revolving nosepiece until it clicks.
- To place the specimen in the field of view, move the stage horizontally back and forth or left and right using the stage movement knobs (7, 8).
- Make sure that the lens does not touch the sample when adjusting the focus: rotate the coarse focusing knob (10) until the specimen is about 3mm (1/8”) away from the lens.
- Turn the fine focusing knob (9) to make the image clear. The fine focusing mechanism allows you to focus on the observed specimen while using high magnifications.
- Change magnification on the screen if necessary. Rotate the diaphragm disk (5) to change the aperture.

Image viewing

KM3 model: To view the microscope image on your smartphone, connect the two devices via Wi-Fi. Use the app to adjust the image and capture photos and videos.

KM5 model: To view the microscope image on your PC, plug one end of the HDMI cable into the corresponding port (4) on the microscope and the other end into your computer. Use the on-screen buttons to adjust the image and capture photos and videos.

KM7 model: Use either of the methods described above.

Main screen menu

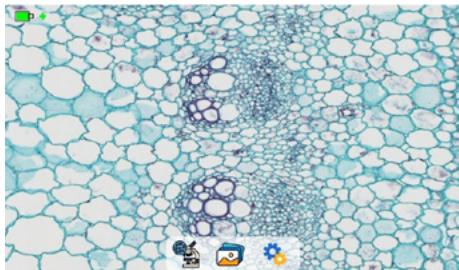
Image	Function
	Battery charge level
	Charging indicator
	Tap to open the microscope menu.
	Tap to open the gallery.
	Tap to open the general settings.
	Tap to view information about all the functions.

Settings

Camera resolution: Tap to select resolution.

Language: Tap to select the interface language (Simplified Chinese, Traditional Chinese, English, Russian, German, French, Japanese, Italian).

USB mode: Tap to select the USB camera mode or Storage mode.

- **USB camera mode:** Connect the USB cable to the port (1) and to your PC. The computer will recognize the microscope as a camera. Open the camera app on your PC to capture photos and videos of the specimen, and save them to your computer.
- **Storage mode:** Connect the USB cable to the port (1) and to your PC. The computer will recognize the microscope as a USB flash drive. Open the folder to view photos and videos saved to the memory card in the microscope, and upload them to your computer.

Storage: Tap to select a USB flash drive or a memory card to store photos and videos.

- **Memory card:** Turn the microscope off and insert the microSD card into the slot at the bottom of the screen until it locks in place with a click. Do not apply excessive force when inserting the microSD card. Try rotating the card if it does not lock in place with ease. Format the card before using it. Turn the microscope on.
- **USB flash drive:** Turn the microscope off and insert the flash drive into the USB port (2). Turn the microscope on.
- If you are using both types of the external storage, tap “Micro SD” or “USB” to switch between them.

Transparency: Use the on-screen slider to adjust the transparency of the menu bar.

Factory oriented: Tap to reset settings.

Microscope menu

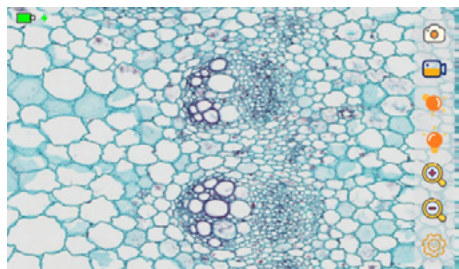
Image	Function
	Tap to take a photo.
	Tap to start recording a video. Tap again to stop recording.
	Tap to turn the upper light on. Use the slider to adjust the brightness. Tap again to turn the upper light off.
	Tap to turn the lower light on. Use the slider to adjust the brightness. Tap again to turn the lower light off.
	Tap to zoom in on the image. Maximum magnification is 3.5x.
	Tap to zoom out on the image.
	Tap to open/close the camera settings menu.

NOTE! The maximum recording time is 20 minutes. Once this time is reached, the current recording will end and a new one will start automatically.

Camera settings menu

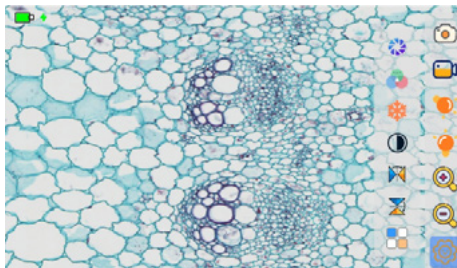
Image	Function
	Tap to turn the auto-exposure function on/off.
	Tap to turn the auto-white balance function on/off.
	Tap to turn the freeze function on/off.
	Tap to turn the grayscale function on/off.
	Tap to flip the image horizontally.
	Tap to flip the image vertically.
	Tap to display or hide a grid and select its color.
	Tap to adjust the exposure and white balance manually.

Gallery menu

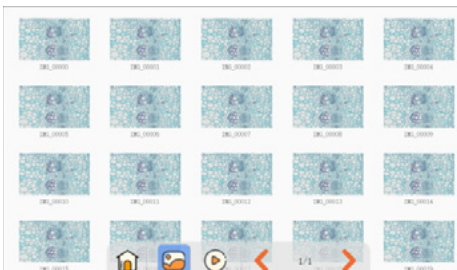
Image	Function
	Tap to exit the preview mode and return to the main page.
	Tap to open the image preview page. Tap the image to enlarge it.
	Tap to open the video preview page. Tap the video to play it.
	Tap to go to the previous page.
	Tap to go to the next page.

Image/video view menu

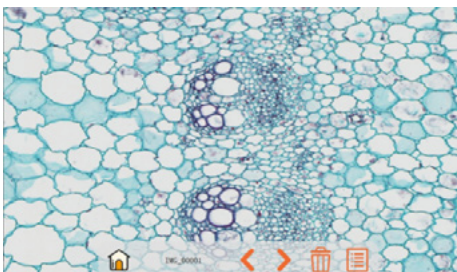
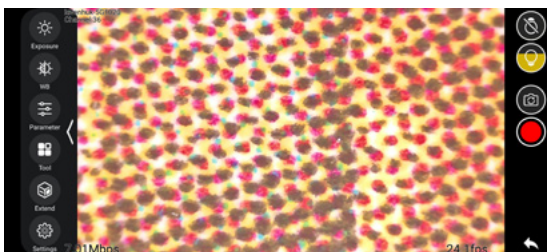
Image	Function
	Tap to exit the single image/video preview mode and return to the main page.
	Tap to go to the previous image/video.
	Tap to go to the next image/video.
	Tap to delete the current image/video.
	Tap to display the current image/video details.

TSViewer app

The TSViewer app provides advanced features:

- Photo/video:** Take photos and videos of a specimen and save it to your device.
- Camera adjustment:** Adjust the exposure, white balance, and resolution.
- Image control:** Use freeze, grayscale, flip, and other functions.
- Illumination:** Control the light sources of the microscope.

Image	Function	Image	Function
	Tap to adjust the exposure.		Tap to set the language, power frequency, and view the software information.
	Tap to adjust the white balance.		Tap to turn the upper light on. Use the slider to adjust the brightness. Tap again to turn the upper light off.
	Tap to adjust the contrast, brightness, etc.		Tap to turn the lower light on. Use the slider to adjust the brightness. Tap again to turn the lower light off.
	Tap to flip, freeze the image, etc.		Tap to switch between photo and video mode.
	Tap to adjust resolution.		Tap to take a photo or record a video. Tap again to stop recording.

S-Viewer software

To work with the photomicroscope on a PC, you need to install the S-Viewer software. The software and the guide are available for download on the official Levenhuk website.

Specifications

	Kelvin Touch KM3	Kelvin Touch KM5	Kelvin Touch KM7
Magnification		53x, 128x, 530x (+ digital zoom 3.5x)	
Microscope head	-	LCD monitor 7" IPS, color 1024x600px, 60fps 2MP camera	LCD monitor 7" IPS, color 1024x600px, 60fps 4MP camera
Photo resolution	2MP	2MP, 5MP	4MP, 8MP
Video resolution	1920x1080px	1920x1080px	1920x1080px; 2560x1440px
Image format		JPEG	
Video format		MP4	
Revolving nosepiece		for 3 objectives	
Objectives, x/aperture		achromatic: 4x/0.10; 10x/0.25; 40x/0.65	
Stage		mechanical stage, stage size: 105x95mm, moving range: 50x20mm	
Condenser		N.A. 0.65, with disc diaphragm	
Focusing mechanism		coaxial coarse & fine focusing knobs	
Light source		LED 0.2W , upper and lower	
Wi-Fi module	5.8GHz	-	5.8GHz
Mobile app	TSViewer	-	TSViewer
App system requirements	iOS 15.6 and above / Android 13 and above	-	iOS 15.6 and above / Android 13 and above
Software	-	S-Viewer	S-Viewer
Software system requirements	-	Windows 7 and above	Windows 7 and above
Storage	-	16GB microSD card (included), up to 512GB	16GB microSD card (included), up to 512GB
Power supply	110-240V via 5V/2A adapter or 1 pc 18650 battery (not included)		

The manufacturer reserves the right to make changes to the product range and specifications without prior notice.

Care and maintenance

Never, under any circumstances, look directly at the Sun, another bright source of light or at a laser through this device, as this may cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS. Take necessary precautions when using the device with children or others who have not read or do not fully understand these instructions. After unpacking your microscope and before using it for the first time, check for the integrity and durability of every component and connection. Do not try to disassemble the device on your own for any reason. For repairs and cleaning of any kind, please contact your local specialized service center. Protect the device from sudden impact and excessive mechanical force. Do not apply excessive pressure when adjusting focus. Do not overtighten the locking screws. Do not touch the optical surfaces with your fingers. To clean the device exterior, use only special Levenhuk cleaning wipes and optics cleaning tools. Do not use any corrosive or acetone-based fluids to clean the optics. Abrasive particles, such as sand, should not be wiped off the lenses, but instead blown off or brushed away with a soft brush.

Do not use the device for lengthy periods of time or leave it unattended in direct sunlight. Keep the device away from water and high humidity environments. Be careful during your observations; always place the dust cover to protect the device from dust and stains when not in use or when being stored. If do not use your microscope for a long time, store the objective lenses and eyepieces separately from the microscope. Store the device in a dry, cool place away from hazardous acids and other chemicals, away from heaters, open fires, and other sources of high temperatures. When using the microscope, try not to use it near flammable materials or substances (benzene, paper, cardboard, plastic, etc.), as the base may become hot during operation and cause a fire. Always disconnect the microscope from a power source before opening the base or changing the bulb. Regardless of the bulb type (halogen or incandescent), give it time to cool down before attempting to replace it, and always replace it with the same type of a bulb. Always use the power supply with the proper voltage, i. e., as specified in the specifications of a microscope. Connecting the instrument to a different power outlet may damage the electric circuitry of the microscope, burn out the bulb, or even cause a short circuit. Children should use the device under adult supervision only. **Seek medical advice immediately if a small part or a battery is swallowed.**

Battery safety instructions

The device is equipped with a rechargeable lithium-ion battery. This avoids frequent battery replacement. Always switch the device off when not in use. If the battery charge is low, please recharge the device in time. Do not overheat the battery. Do not discharge the battery completely. Keep batteries out of the reach of children, to avoid risk of ingestion, suffocation, or poisoning. Utilize used batteries as prescribed by your country's laws.

Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a **lifetime** warranty against defects in materials and workmanship. Lifetime warranty is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for **six months** from date of retail purchase. The warranty entitles you to free repair or replacement of the Levenhuk product in any country where a Levenhuk office is located if all warranty conditions are met.

For further details please visit our web site: levenhuk.com/warranty

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch.

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Levenhuk® is registered trademark of Levenhuk, Inc.
© 2006-2026 Levenhuk, Inc. All rights reserved.
Levenhuk Inc. (USA): 6021 Catlin Dr., Tampa, FL 33612, USA,