# Levenhuk Kelvin TMC20 LED Microscope for Smartphone

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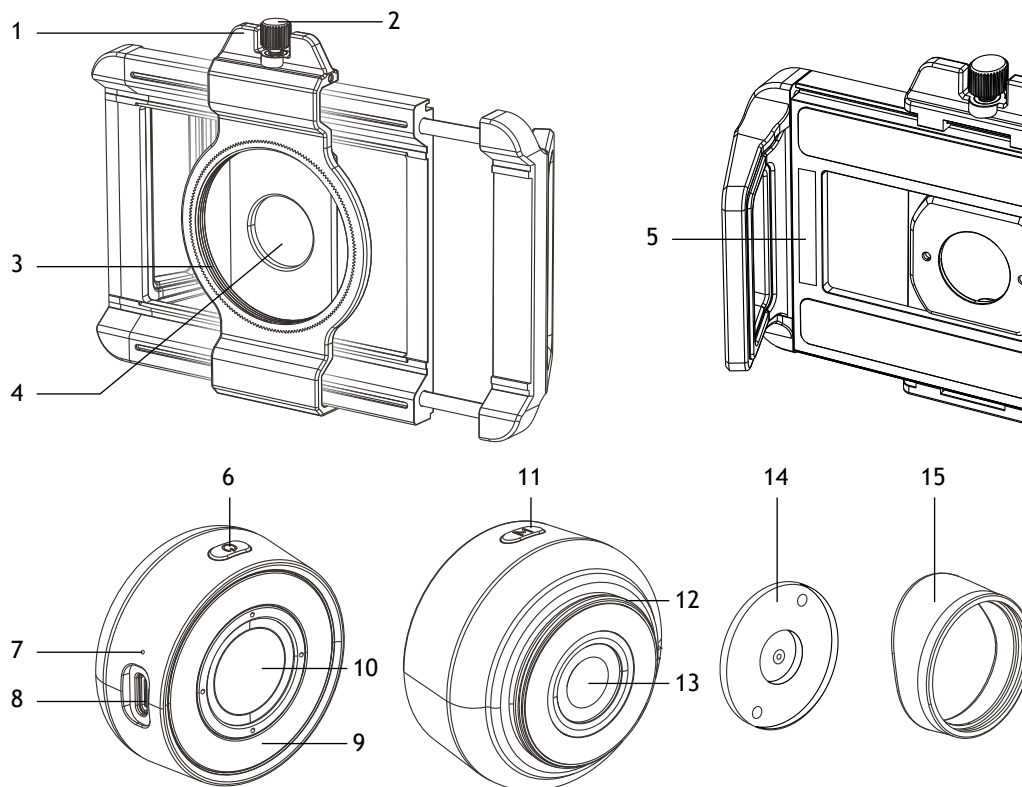
# EN Levenhuk Kelvin TMC20 LED Microscope for Smartphone

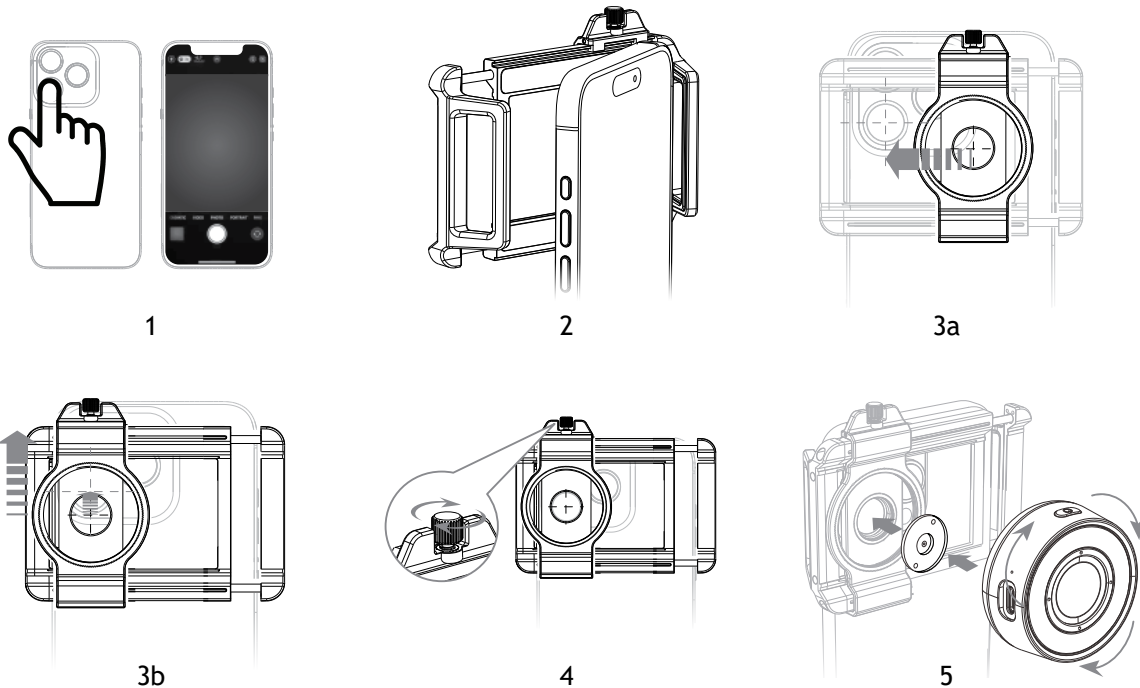
**Caution! Never look directly at the Sun through this device, as this may cause permanent eye damage and even blindness.**

The kit includes: LED microscope, phone-to-lens mounting bracket, USB Type-C cable, eyecup, field of view diaphragm, EVA pads (9 pcs), pouch, cleaning cloth, user manual, and warranty.

## Charging the device

Connect the USB Type-C cable to the USB Type-C slot (8) and power bank, PC, or power adapter (not included). The power indicator (7) will light red. Charging takes approximately 90 minutes. When charging is complete, the power indicator will light green.

## Getting started

- Find the main camera on your smartphone. To do so, open the camera app on your smartphone and cover each lens on the back one by one with your finger. The module that makes the image on the screen go black when covered is the main camera (Fig. 1).
- Gently push your smartphone into the phone-to-lens mounting bracket (Fig. 2). To precisely adjust the mounting bracket to the thickness and width of your phone, use the soft EVA pads (included).
- Move the lens mount (1) horizontally and vertically to align the lens hole (4) with your smartphone's main camera (Fig. 3a, 3b).
- Tighten the fixing screw (2) (Fig. 4).
- Install the field diaphragm (14) directly into the threaded interface (3), and then screw the threaded interface (12) clockwise (Fig. 5).

**! The field diaphragm is not fixed. Take care not to drop it when installing and removing.**

You may now use the device to observe objects, take photos, and record videos using your smartphone.

## Usage

- Press and hold (6) to turn the device on. The LED illumination (9) will light up.
- Press (6) to change the brightness level (low/middle/high).
- Press (11) to change the LED illumination mode. You may choose from 2 color temperatures (cool or warm light) in 3 illumination ranges (full circle / left half circle / right half circle) in each of 3 brightness modes (low/middle/high), or you may select the UV full circle mode.
- Press and hold (6) to turn the device off.

**! If there are uniform dark corners around the image, adjust the digital zoom on your smartphone. If the dark corners look uneven, check that the lens hole (4) is aligned with your smartphone's main camera before trying again.**

You may also observe the objects only using the microscope with the eyecup (15) without installing the device on your smartphone. Please note that this way, taking photos or recording videos is not available.

## Specifications

Magnification	120x
Focusing distance	10mm
Objective lens diameter	22mm
Optics material	optical glass
Optics coating	fully multi-coated
Brightness level	3 modes
LED illumination	19 modes
Compatibility with smartphone models	iPhone, Android; smartphone width: 70–100mm (2.8–3.9"), smartphone thickness: ≤ 20mm (0.8")
EVA pads	24x4.5x2mm (3 pcs) 24x4.5x3mm (3 pcs) 24x4.5x4mm (3 pcs)
Operating temperature range	0... +45° C (+32... 113° F)
Power supply	3.7V, 110mA·h rechargeable Li-ion battery; 5V, 1A DC adapter (not included); USB Type-C cable (included)
Battery duration	120 min

The manufacturer reserves the right to make changes to the product range and specifications without prior notice.

## Care and maintenance

Take the necessary precautions when using the device with children or others who have not read or who do not fully understand these instructions. Do not try to disassemble the device on your own for any reason. Protect the device from water, dust, sudden impact and excessive mechanical force. Do not touch the optical surfaces with your fingers. Clean the lens surface with compressed air or a soft lens cleaning cloth. Never attempt to operate a damaged device or a device with damaged electrical parts! Store the device in a dry cool place. **If a part of the device or battery is swallowed, seek medical attention immediately.**

## Battery safety instructions

The device is equipped with a rechargeable lithium-ion battery. This avoids frequent battery replacement. Always switch the device off when not in use. If the battery charge is low, please recharge the device in time. Do not overheat the battery. Do not discharge the battery completely. Keep batteries out of the reach of children, to avoid risk of ingestion, suffocation, or poisoning. Utilize used batteries as prescribed by your country's laws.

## Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars, and other optical products, except for their accessories, carry a **lifetime warranty** against defects in materials and workmanship. A lifetime warranty is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for **six months** from the purchase date. The warranty entitles you to the free repair or replacement of the Levenhuk product in any country where a Levenhuk office is located if all the warranty conditions are met.

For further details, please visit: [levenhuk.com/warranty](https://www.levenhuk.com/warranty)

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch.

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Levenhuk® is registered trademark of Levenhuk, Inc.
© 2006–2026 Levenhuk, Inc. All rights reserved.
Levenhuk Inc. (USA): 6021 Catlin Dr., Tampa, FL 33612, USA,